

Product Safety Information

Safety cabinet type 90 Asecos Q-CLASSIC-90 asecos, width 1200 mm, variant 1

Manufacturer/Responsible Person: asecos GmbH, Weiherfeldsiedlung 16-18, 63584 Gründau, Deutschland

E-Mail: info@asecos.com

Please note the technical data of the products!

Article Numbers: 132011, 132010, 132009, 132008, 132007, 132006, 132005

Safety Instructions:

- Regular maintenance and functional testing of the safety storage cabinet by a competent person is required (at least once a year).
- The cabinet is designed exclusively for the storage of hazardous substances in accordance with TRGS 510. Read the manufacturer's safety instructions carefully and observe them.
- In the event of fire, the door closes automatically due to the integrated thermal release. Additional fire protection measures must be taken into account.
- Do not block the exhaust air connection and maintain safety distances to avoid hazardous vapours.
- Place the safety storage cabinet on stable surfaces and level out uneven floors to ensure stability.
- Maximum load: shelf - 75 kg, pull-out tray - 60 kg; do not exceed.
- Do not overload the doors to prevent the locking mechanism from failing.
- Secure the cabinet during transport to prevent injury or damage.
- Take particular care when storing flammable substances.
- Observe the safety and warning notices and regulations in accordance with TRGS 510 on the product and in the operating instructions.
- Storage temperature should comply with the specifications for hazardous substances.
- Use safety containers such as floor drip pans to prevent leakage.