

Product Safety Information

Safety cabinet type 90 asecos, W 1196 x D 616 x H 1968 mm, 3 shelves & drip tray asecos, l.grey RAL 7035 or s.yellow RAL 1004

Manufacturer/Responsible Person: asecos GmbH, Weiherfeldsiedlung 16-18, 63584 Gründau, Deutschland

E-Mail: info@asecos.com

Please note the technical data of the products!

Article Numbers: 22099, 22098

Safety Instructions:

- Ensure that the cabinet is properly and securely locked to prevent access to hazardous substances. Use the profile cylinder lock to prevent unauthorised access.
- In the event of fire, the doors close automatically. Check the door hold-open system, which has a self-closing function in the event of a fire.
- Observe safety measures when handling flammable liquids and comply with the relevant regulations in accordance with TRGS 510.
- The cabinet must be placed on a stable, level surface to prevent it from tipping over.
- Ensure that the storage levels are loaded in accordance with the permissible loads (tray shape: max. 75 kg, base tray: max. 150 kg).
- Check the cabinet regularly for damage or wear and tear and repair or replace if necessary.
- Fire hazard: The cabinet is intended for the storage of flammable liquids. Therefore, comply with all relevant safety standards for the use of such a cabinet.
- Safety ventilation: Check that the exhaust air connection is working properly.
- Warnings for safe use: Keep the operating doors wide open during use and ensure that the cabinet is not overloaded.