

Product Safety Information

Chemical cabinet asecos CS-CLASSIC, hinged doors, W 810 x D 520 x H 1105 mm

Manufacturer/Responsible Person: asecos GmbH, Weiherfeldsiedlung 16-18, 63584 Gründau, Deutschland

E-Mail: info@asecos.com

Please note the technical data of the products!

Article Numbers: 463112, 463111, 463110, 463109, 463108, 463107, 463106, 463105, 463104, 463103, 463102, 463101, 463100, 463099

Safety Instructions:

- Use suitable personal protective equipment such as gloves and safety goggles for the safe handling of hazardous substances.
- Ensure that the chemical cabinet is located in a well-ventilated area and that the ventilation slots are not blocked. Ensure proper installation and regular maintenance of the ventilation system to safely dissipate hazardous vapours.
- Use a technical exhaust air system to minimise the risk of explosion.
- The chemical cabinet may only be used by trained and authorised personnel to prevent unauthorised access. Use the integrated cylinder lock for this purpose.
- Keep the cabinet locked when it is not in use to prevent unauthorised access.
- Regularly check the tightness of the base drip tray and the entire cabinet for leaks and other damage. In the event of damage, take immediate action or contact the manufacturer.
- Ensure that the maximum permissible load capacity of the shelves (60 kg) is not exceeded.
- Store and dispose of flammable materials and hazardous chemicals safely and in accordance with local regulations.
- Only suitable for indoor use.
- Read and follow all safety information, maintenance manuals and instructions for use, which should be provided in an easily accessible and understandable language.
- The chemical cabinet should only be used for suitable chemicals and not as storage space for other items.